



### PRODUCT DESCRIPTION

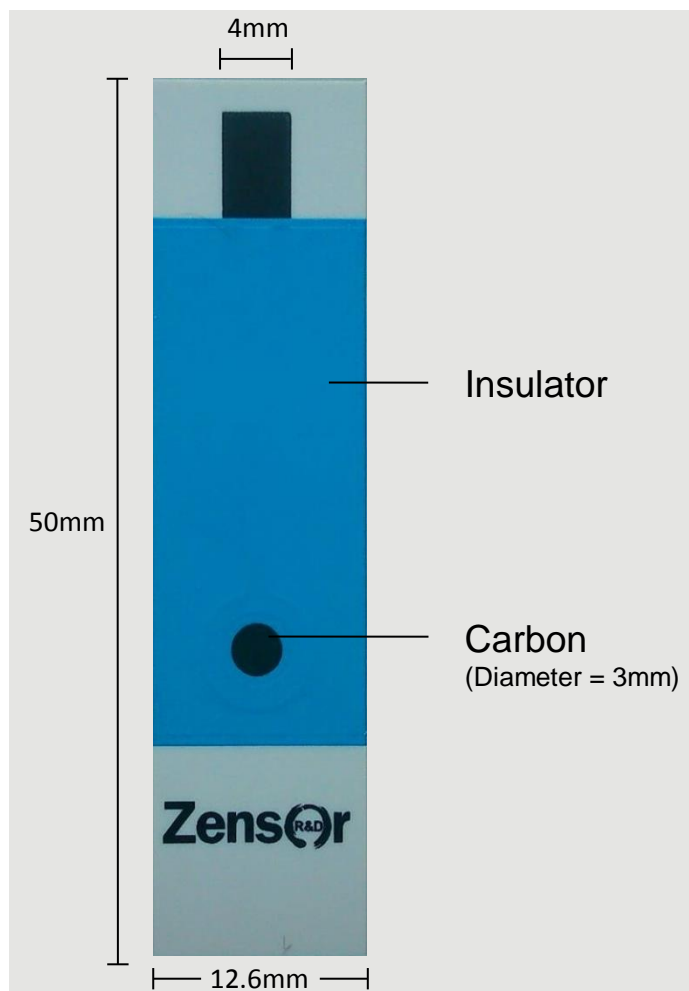

### Before Use (Recommend)

- A** You can wash DI water with several times.

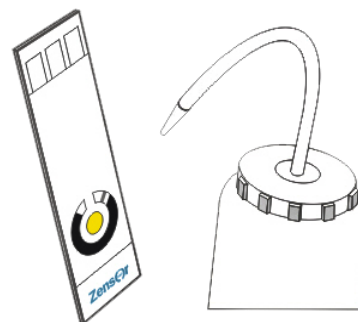

- B** Or put electrode of SPE into DI water or buffer solution of you prepared at 5~10 min.

- C** Or put electrode of SPE in 0.1M phosphate buffer solution (pH7.0) and then operation of cyclic voltammetry (CV) with sever segments.

### DESCRIPTION FOR USE

1. SE101 is supplied ready for use under acid ( $< 2 \text{ M H}_2\text{SO}_4$ ) and basic media ( $< 1 \text{ M NaOH}$ ).
2. Avoid bending hardly as operating.
3. Should Keeping apart from organic media.

### Storage

Store product in a dry location apart from sunlight at optimal temperature of 20 to 40 °C.